

Discovery MX10 Professional Incubator

The Discovery MX10 is designed to incubate up to 8 (10 screw top type) Dip slides and can be powered from the mains via the adaptor supplied or via a cigar lighter cable in a vehicle.

For mains use select the applicable plug for your location and clip in the adaptor, Power input is located on the LH side of the Discovery case, to use in vehicle is simply a matter of changing cable.

A QUICKSET guide point is shown on the temperature control at 37°C, these should be used for reference, each quarter turn of the temperature control equates to 7°C+/- from this point so for example roughly midway (12.00) would be around 30°C – ideal for L8 dip slide applications.

The actual temperature must always be taken against the thermometer inside the unit as with all incubators large variants in ambient temperatures can influence internals slightly
Prior to loading samples it is recommended that the incubator should be warmed up for a period of 30 minutes to stabilize on mains power

The mains power pack features automatic overload protection and the 12v cigar lighter cable has a 1.5a QB fuse in the large plug end. Please refer any service or repair to Lynd Products Ltd or your supplier.

As with all our Incubator products this Discovery MX10 benefits from a 24 month warranty against faulty workmanship, date of invoice is deemed date of supply.