

## Calcium Hardness TC

UFI: QVGE-ER1N-X413-4KD9



### Danger

**H302** - Harmful if swallowed

**H314** - Causes severe skin burns and eye damage

**H332** - Harmful if inhaled

**P260** - Do not breath dust/fume/gas/mist/vapours/spray

**P280** - Wear protective gloves/protective clothing and eye eye/face protection

**P303 + P361 + P353** - IF ON SKIN (or hair): Take off immediatly all contaminated clothing. Rinse skin with water [or shower]

**P305 + P351 + P338** - IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing

**P310** - Immediately call a POISON CENTER or doctor

# Test Kit

CaCO<sub>3</sub> (10 - 500 mg/l)



## Calcium Hardness TC

**NOTRUF/EMERGENCY: +49 (0)89-19240 (24h)**

Water-i.d. GmbH  
Daimlerstr. 20  
D-76344 Eggenstein  
Tel. +49 (0) 721 - 782029-0  
Fax +49 (0) 721 - 782029-11  
[www.water-id.com](http://www.water-id.com)  
[info@water-id.com](mailto:info@water-id.com)



**Calcium Hardness** Tablet-count method



**Calciumhärte** Tabletten-Zählverfahren



**Durété calcique** Methode de comptage à pilule



**Durezza del calcio** Metodo di conteggio de pastiglie









**Dureza cálcia** Método Contador de tabletas




**Calciumhårdhed** Tablettællings-metoden



**Calcium hardheid** Tablet-telmethode

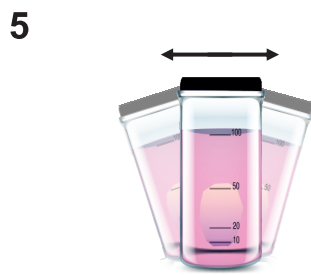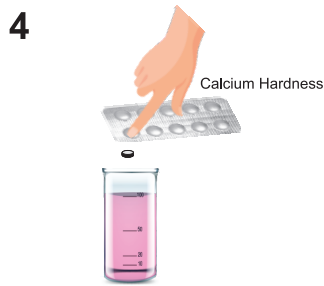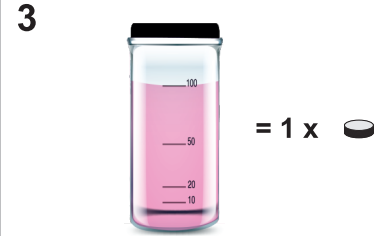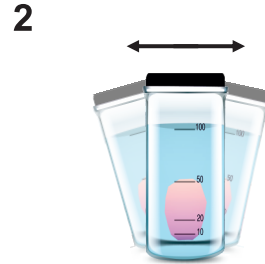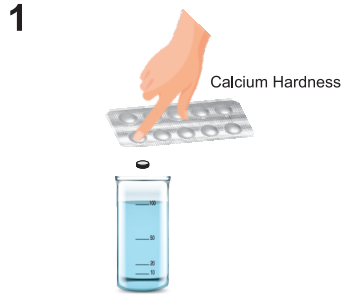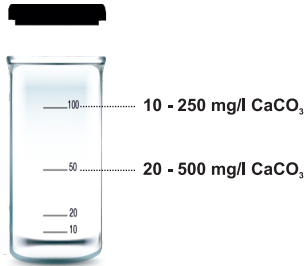

50 ml: (  x 40) - 20 = mg/l CaCO<sub>3</sub>

100 ml: (  x 20) - 10 = mg/l CaCO<sub>3</sub>

	mmol/l	mval/l	°dH	°e	°f
1 mg/l CaCO <sub>3</sub>	0.01	0.02	0.056	0.07	0.1

Alkaline Earth Ions  
Erdalkali-Ionen  
Ioni alcalino-terreux  
Ioni alcalino-terrosi  
Iones alcalina terrosos  
Löes alcalino-terrosos  
Ioner fra alkalisk jord  
aardalkali-ioner

